




## Battery activation


Open the top cover to access the cooling unit.

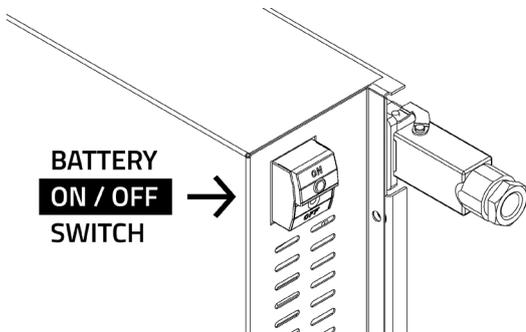

Turn on the On the battery switch.

## Dialog connection

Plug the device to the mains and turn it on.

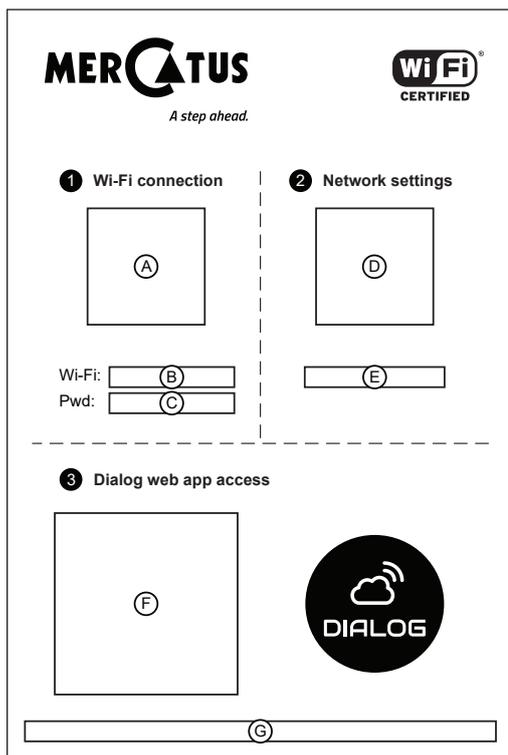

For smartphones and tables:

1. Scan QR code (A)
2. Connect Wi-Fi (B) and enter PWD (C)
3. Scan QR code (D) to access network settings
4. Insert your network SSID and PWD
5. Go to Wi-Fi and connect it to your local Wi-Fi
6. Scan QR code (E) and register on Dialog web app
7. Now onwards your device is connected, sign in Dialog to access all data

For laptops and desktops:

1. Turn on the Wi-Fi connect it to (B) and enter PWD (C)
2. Open browser and access (E)
3. Insert your network SSID and PWD
4. Go to Wi-Fi and connect it to your local Wi-Fi
5. Access (G) and register on Dialog web app
6. Now onwards your device is connected, sign in Dialog to access all data

For more information, see Dialog sync Manual.

## Volt-free contact connection

The connection allows to transmit the following alarms: High temperature; Low temperature; Open door; Power failure. To consult the connection diagram, see User Manual.

## Antifreeze module management

This is a safety feature that prevents temperature from dropping below the set minimum. Acura models feature a digital protection while Exige models feature a mechanical thermostat. Factory setting is at 2,5°C, to edit this setting see Controller Manual.

## Useful information

For more information, see the instruction manual:  
<https://bioscience.mercatus.pt/en/resources>