
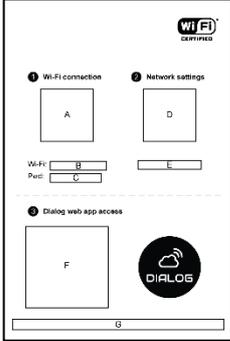
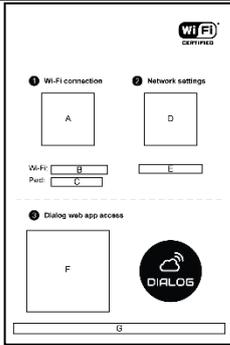



Step-by-Step connection to Dialog IoT

- 1 Connect the plug to the socket to turn on the equipment. In the first connection, you must be physically close to the equipment.


- 2  In the top left corner of the device's interior you can find the label for sync. Alternatively, you can consult the additional label that is sent along with the equipment documents.

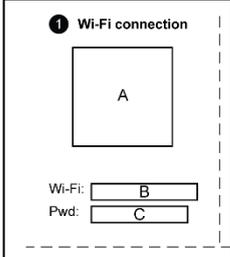



- 3  Scan QR code **A** to connect to Wi-Fi. Connect to the Wi-Fi network **B** with the name and serial number of your device. Ex: Dialog Sense <Model> <serial number>. Enter the provided password **C**.

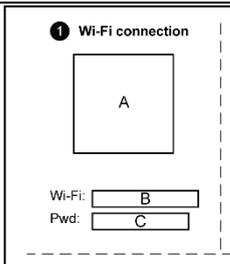


- 4  If a similar message appears, press confirm or connect. Wait a few moments and proceed when the Wi-Fi **B** is connected.

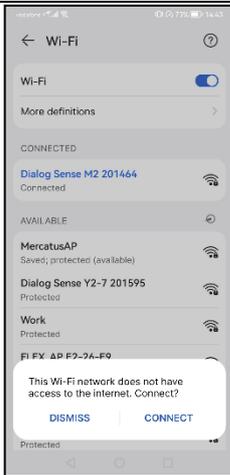

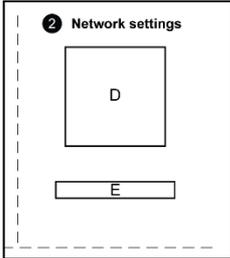- 5  Scan the QR code **D** to connect to the network or enter **http://192.168.2.1** in your browser.

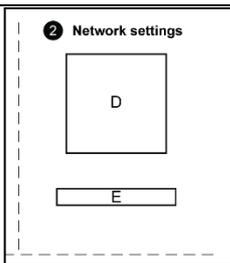


6

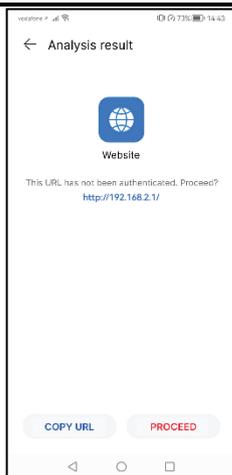

If a similar message appears, press confirm or proceed.

7

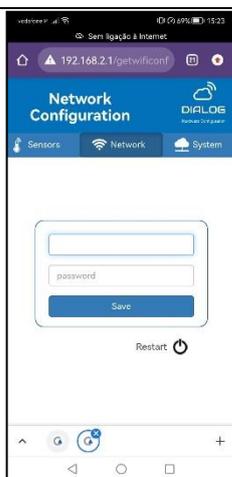

You will be taken to this page, go to **Network** tab. Here you must enter your network data: SSID * and password. After entering the data, press the **Save** button.

* The SSID is the name of your Wi-Fi network.


8



On the same page, select the **System** tab.


Wait a few moments until data starts to appear. After that, your equipment is connected to the network, you can close the page.

9

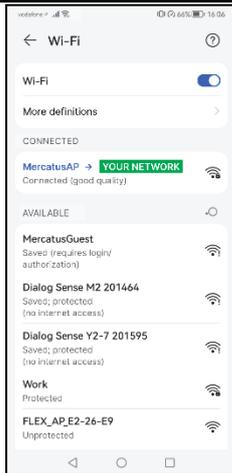

Go to your Wi-Fi settings and connect to your local Wi-Fi network.

10

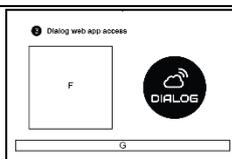


Scan the QR code **F** to connect to the network or write the link **G** in your browser.

11




If a similar message appears, press confirm or proceed.

12

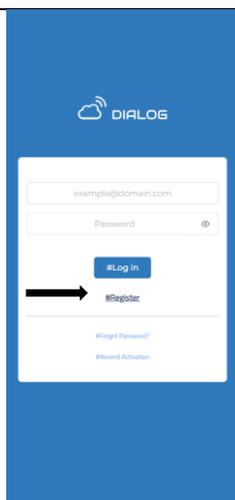

You will be directed to the next page.
Select **Register** to create your Dialog account.

13

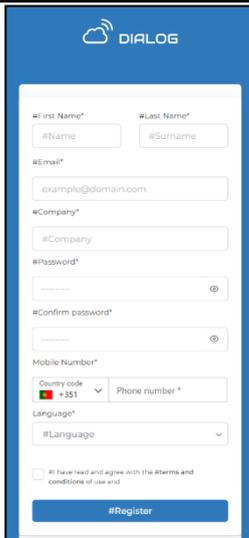A screenshot of a mobile registration form for Dialog. The form is titled 'DIALOG' and contains several input fields: '#First Name*' and '#Last Name*' (with sub-fields '#Name' and '#Surname'), '#Email*' (with the example 'example@domain.com'), '#Company*', '#Password*' (with a strength indicator), and '#Confirm password*'. Below these is the 'Mobile Number*' section with a 'Country code' dropdown (set to '+351') and a 'Phone number *' field. There is also a 'Language*' dropdown menu. At the bottom, there is a checkbox for '#I have read and agree with the terms and conditions of use and' and a blue '#Register' button.

Fill in the registration details. After completing the form, click on the **Register** button to proceed. Introduce a payment method and enjoy the free trial.

14

Go to your email inbox. Have received an email from noreply.dialog@mercatus.pt. Click on the provided link to active and login to your Dialog account.

15

From now on, you can access Dialog by simply entering your login credentials from any device with a Wi-Fi connection.

To access Dialog use the link iot.mercatus.pt.